




## Product Codes

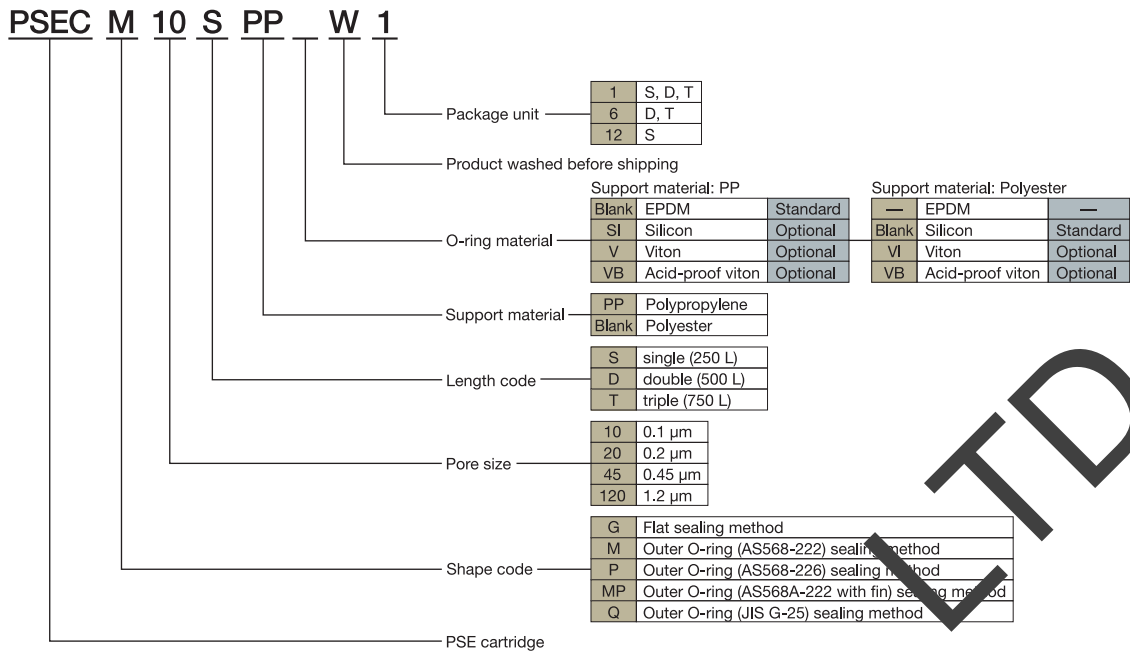

## Sterilization Process

The AstroPore PSE Cartridge W can be sterilized repeatedly. Sterilization at specified intervals is required for maintaining reliable and consistent bacterial removal and filtration performance.

Hot water sterilization : 90°C


Chemical sterilization :

- Hydrogen peroxidized 5%
- Formalin 1%
- Sodium hypochlorite 10ppm (Chlorine water)

Steam sterilization : 121°C \* 30mins \* 30 cycles  
(For more cycles requirements, please contact sales representative)

## Types

The following three types are available depending on the form of sealing, G, M, P and MP types.

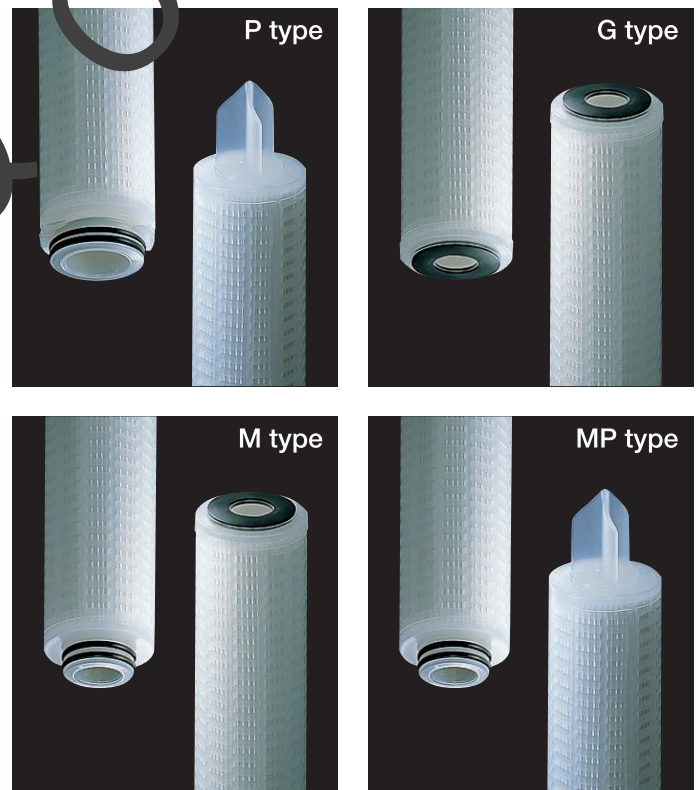

**FUJIFILM**

FUJIFILM Corporation

<http://www.fujifilm.com/products/microfilter/>

TEL: +81-3-6271-3008

誠風企業股份有限公司

DEOVOS & Co Ltd

10485台北市建國北路2段66號3樓之3

T.02-2516-2881 F.02-2516-2883

[www.devos.com.tw](http://www.devos.com.tw)